

QUANTOFIX® Ascorbic acid

Pack content:

1 aluminum container with 100 test strips

Measuring range:

Visually	Reflectometrically
50–2000 mg/L	25–1000 mg/L

Color gradation:

0 · 50 · 100 · 200 · 300 · 500 · 700 · 1000 · 2000 mg/L

Hazard warnings:

This test does not contain hazardous substances that must be labelled.

General indications:

Remove only as many test strips as are required. Close the container immediately after removing a strip. Do not touch the test field.

Instructions for use:

1. Dip the test strip into the tests solution for 10 s.
2. Shake off excess liquid.
3. Wait 30 s.
4. Compare test field with the color scale.

If ascorbic acid is present, the test field turns green to blue.





Disposal:

Information regarding disposal can be found in the safety data sheet. You can download the SDS from www.mn-net.com/SDS.

Storage:

Avoid exposing the strips to sunlight and moisture. Keep container cool and dry (storage temperature 4–30 °C).

If correctly stored, the test strips may be used until the use-by-date printed on the packaging.

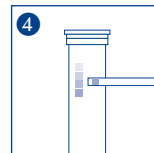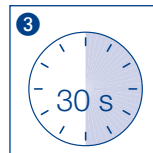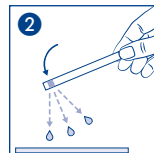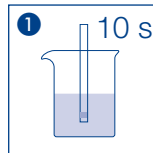

CTL SCIENTIFIC SUPPLY CORP. 1016-3 Grand Boulevard, Deer Park, NY 11729

Tel: 631-242-4249

Web: www.ctlscientific.com

Manufacturer: Macherey-Nagel GmbH & CO. KG

Rev: 2025-03