

# QUANTOFIX® Ascorbic acid

Pack content:

1 aluminum container with 100 test strips

Measuring range:

Visually  
50–2000 mg/L

Reflectometrically  
50–1000 mg/L

Color gradation:

0 · 50 · 100 · 200 · 300 · 500 · 700 · 1000 · 2000 mg/L

General indications:

Remove only as many test strips as are required. Close the container immediately after removing a strip. Do not touch the test field.

Instructions for use:

1. Dip the test strip into the tests solution for 10 s.
2. Shake off excess liquid.
3. Wait 30 s.
4. Compare test field with the color scale.

If ascorbic acid is present, the test field turns green to blue.


Disposal:

Used test strips can be placed in the normal household waste.

Storage:

Avoid exposing the strips to sunlight and moisture. Keep container cool and dry (storage temperature not above +30 °C).

If correctly stored, the test strips may be used until the use-by-date printed on the packaging.

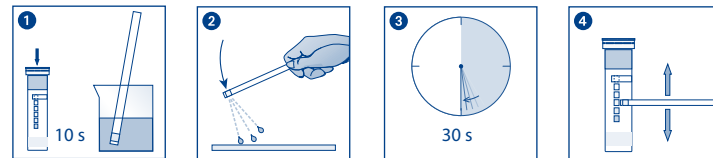

## Product data and ordering information

Type	Semi-quantitative test strips
Presentation	Tube of 100 test strips
Color reaction	yellow to green-blue
Gradation	0 · 50 · 100 · 200 · 300 · 500 · 1000 · 2000 mg/l (ppm) Ascorbic acid / Vitamin C
REF	91314