

# pH-Fix 0 – 14

## Test strips for measuring pH values in the range of 0 – 14



Please refer to instructions for use

**REF** 92110

### Pack content:

100 test strips

1 reference color scale pH 0 – 14

### Instructions for use:

1. Dip test strip for 1 s with all test fields in the sample solution.
2. Shake off excess sample solution.
3. Compare with color scale and read off the corresponding pH-value.



### Note:

If sample solution is weakly buffered leave test strip in solution until color no longer changes.

### Storage:

Avoid exposing the test strips to sunlight and moisture. Store the container in a dry place (storage temperature 4 – 30 °C).

### Disposal:

Used test strips may be disposed of as domestic waste.



Please read instructions for use!

**REF**

Item number

CTL SCIENTIFIC SUPPLY CORP. 1016-3 Grand Boulevard, Deer Park, NY 11729

Tel: 631-242-4249

Web: [www.ctlscientific.com](http://www.ctlscientific.com)

Manufacturer: Macherey-Nagel GmbH & CO. KG

Rev: 2025-03