


pH-Fix 4.5–10.0

Test strips for measuring the pH values in the range of pH 4.5–10.0

 Please refer to instructions for use

REF 92120

Pack content:

100 test strips

1 reference color scale pH 4.5–10.0

Application:

The pH-Fix test strips pH 4.5–10.0 are semi quantitative indicator test strips for many applications.

For quality control use commercially available pH buffer solution.

Instruction:

1. Dip the test strip with all its fields into the sample solution for 1 sec.
2. Shake off excess liquid.
3. Compare the colors with the color chart and read off the pH value which matches closest.



Remark:

In weakly buffered solutions, keep the test strip immersed for a longer period of time until color consistency of all test fields.

Storage:

Avoid exposing the strips to sunlight and/or moisture. Keep container cool and dry (store between 4–30 °C).

Disposal:

Used test strips may be disposed of as domestic waste.

 Please read instructions for use!

 Permitted storage temperature range

 Use by

LOT Batch identification

REF Item number

 Manufacturer

 Do not reuse