

pH-Fix 6.0 – 10.0

Test strips for measuring pH values in the range of 6.0 – 10.0



Please refer to instructions for use

REF

92122


Pack content:

100 test strips


1 reference color scale pH 6.0 – 10.0

Instructions for use:


1. Dip test strip for 1 s with all test fields in the sample solution.
2. Shake off excess sample solution.
3. Compare with color scale and read off the corresponding pH-value.

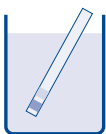

1

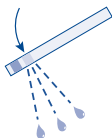

2

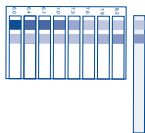

3

Note:

If sample solution is weakly buffered leave test strip in solution until color no longer changes.

Storage:

Avoid exposing the test strips to sunlight and moisture. Store the container in a dry place (storage temperature 4 – 30 °C).

Disposal:

Used test strips may be disposed of as domestic waste.



Please read instructions for use!

REF

Item number

CTL SCIENTIFIC SUPPLY CORP. 1016-3 Grand Boulevard, Deer Park, NY 11729

Tel: 631-242-4249

Web: www.ctlscientific.com

Manufacturer: Macherey-Nagel GmbH & CO. KG

Rev: 2025-03