


# pH-Fix 3.6 – 6.1

## Test strips for measuring the pH values in the range of pH 3.6 – 6.1

 Please refer to instructions for use

**REF** 92130

**Pack content:**

100 test strips

1 reference color scale pH 3.6 – 6.1

**Instruction:**

1. Dip the test strip with all its fields into the sample solution for 1 sec.
2. Shake off excess liquid.
3. Compare the colors with the color chart and read off the pH value which matches closest.



**Remark:**

In weakly buffered solutions, keep the test strip immersed for a longer period of time until color consistency of all test fields.

**Storage:**

Avoid exposing the strips to sunlight and/or moisture. Keep container cool and dry (store between 4 – 30 °C).

**Disposal:**

Used test strips may be disposed of as domestic waste.

 Please read instructions for use!

**REF** Item number