

Chlorine Test

Chlorine Test is a test paper for the semi-quantitative determination of chlorine in the range of 10-200 ppm.

Color reaction:

In the presence of chlorine the white test paper changes color to blue-violet.

Instructions for use:

Dip test strip **briefly** into the test solution and shake off excess of liquid. Compare reaction color **immediately** with color scale.

Remark:

If the test paper is discolored after dipping, the chlorine concentration is higher than 200 ppm. In this case the result should be ignored.

Storage:

Avoid exposing the test paper to sunlight and moisture. Store the container below +30°C in a dry place.

Product data and ordering information

Type	special test paper
Presentation	Reel of 5 m length
Color reaction	white to dark-blue
Range	0 · 10 · 50 · 100 · 200 mg/l (ppm) Cl ₂
REF	90709